


Components list

Thank you for choosing our products. Please read the assembly steps and user manual carefully before use.

If you have any problems with the use and assembly, please feel free to contact us.

We will reply within 24H (there may be a delay on holidays). We appreciate the opportunity to serve you, your satisfaction is our utmost priority!

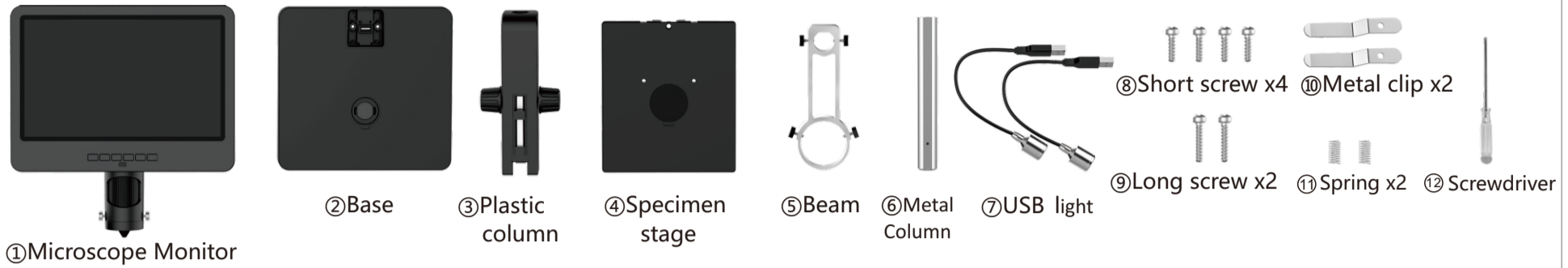

Assembly steps

Step1

Insert ③ into ②.
*Direction as shown

Step2

Fix short screws ⑧ into the base ②.

Step3

Insert ④ into ③. Fix a short screw ⑧ into the specimen stage ④.

Step4

Insert ⑥ into ③. Assemble as shown in the figure. Fix a short screw ⑧.

Step5

Slot ⑤ to ⑥. Tighten the screws.

Step6

Unscrew the screws and remove the lens.
*Be careful not to drop the lens.

Step7

Insert the Protrusion into the slot. Assemble the lens and tighten the screws.

Step8

This step is only required when using lens A to observe coins, insects, stones, etc., and it is not required when using the Lens B&C to observe slides.

Assemble the USB light ⑦.

Option

This step is only required when observing live insects such as butterflies, and it is not required for other purposes.

Use long screws ⑨ for metal clips assembly.